

Applied Sciences

→ UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.



Applied Sciences

UV & Sunblock






Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Objective

The purpose of this activity is to study the level of ultraviolet radiation resulting from the intervention of a sunlight beam through different types of filters, creating a hypothesis and proceeding to test it using the Labdisc UV sensor.

The sun has always played a unique role in our environment, giving energy to living beings and directly warming up the planet. Human beings have always had to coexist with benefits and damages emerging from the presence of the sun. However, today we have more knowledge about the damaging effects of too much exposure to the sun.

UV index table

UV-index Level	Exposure Category	Color Code	Precautions
2 or less	Low		You can safely be outdoors with minimal protection.
3 or 5	Moderate		Wear a hat, sunscreen, sunglasses or seek shady areas.
6 to 7	High		Wear a hat, sunscreen, sunglasses or seek shady areas. Stay indoors between 10am and 2 pm (11-3 daylight saving time).
8 to 10	Very High		Stay indoors as much as possible. otherwise use all precautions above.
11 or higher	Extreme		As above

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Introduction and theory

?

How could you physically define the sun radiation?

?

Describe the positive and negative effects produced by the sun on Earth. Do you know methods to prevent negative effects from sun exposure?

Carry out the experiment activity with your class so that at the end you'll be able to answer the following question:



?

What is the precise role of skin products and eyewear that protect against the sun?

Measuring and comparing the level of ultraviolet radiation, through different types of filters.


Theoretical

The sun radiation is electromagnetic (EM) energy, which is transmitted through a vacuum. The EM radiation is made up of packets of energy or photons, which travel from the sun in a wave-like pattern. There is a scientific scale where the EM radiation is classified according to the wavelength, frequency and energy of the radiation: The EM spectrum. The UV radiation is a form of energy placed between the visible spectrum and X-rays. It is known that UV rays help the body to produce vitamin D3 and are also very important to create oxygen through photosynthesis. However prolonged exposure could cause eye and skin damage as a short-term consequence and skin cancer as a side effect.

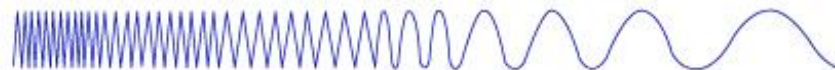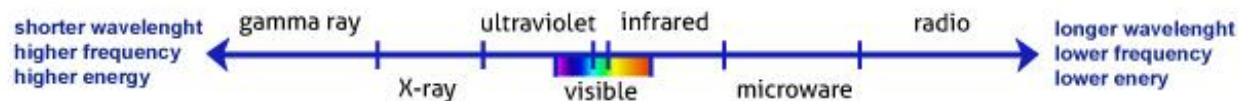

Sunlight brings three types of UV rays: UVA (longest wavelength), UVB and UVC (shortest wavelength). The UV radiation that touches us is made up of UVA and UVB rays, whereas UVC rays are absorbed by the ozone layer. Part of the UVB rays are also rejected by this layer. While the UVA reaches a deeper skin layer. UVB is the most harmful radiation to the eye.

The sun protectors absorb and reflect the UV radiation range and other electromagnetic waves, acting as a filters. Sunglasses and sunscreens have chemicals that perform this function, thus Preventing direct passage of the UV energy to the eyes or skin.

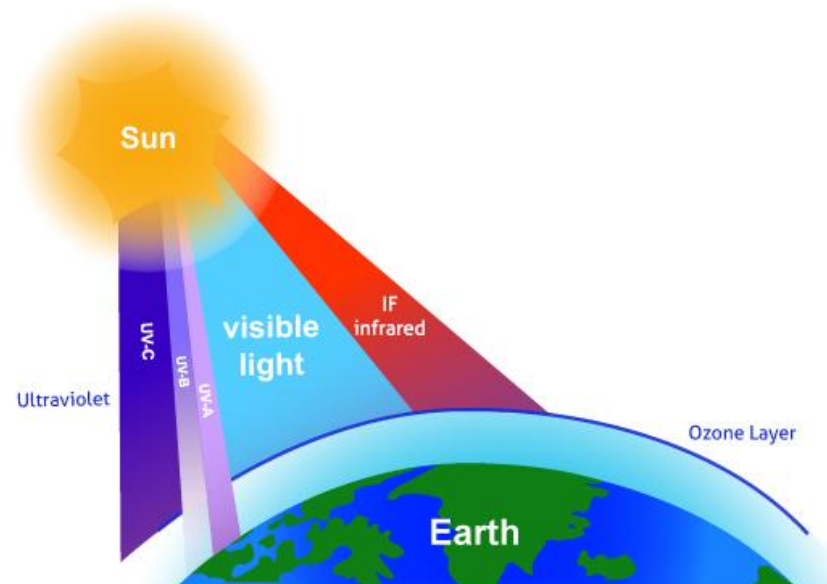

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Introduction and theory

Now students are encouraged to raise a hypothesis which must be tested with an experiment.



How effective are the classical methods to prevent sun damage? Are there differences between these types of sun filters?

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Activity description

Students will measure ultraviolet radiation from the sun, testing different types of sun filters, such as sunblocks (SPF 15 and 30) and sunglasses. They will evaluate these objects and substances according to the experimental results, using the Labdisc UV sensor.

- 1 Labdisc Enviro
- 2 1 plastic transparent film (10 x 6 cm)
- 3 1 cardboard (10 x 6 cm)
- 4 Liquid silicone
- 5 2 pairs of sunglasses
- 6 2 sunblocks (FPS 15 and 30 preferably)

Applied Sciences

UV & Sunblock


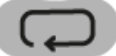

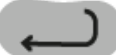

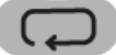






Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Resources and materials



a. Labdisc configuration

To collect measurements with the UV sensor, the Labdisc must be configured following these steps:

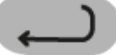- 1 Turn on the Labdisc pressing  .
- 2 Press  and select "SETUP" by pressing  .
- 3 Now select the option "SET SENSORS" with  , and choose "UV". Then press  .
- 4 Once you have done that, you will be back at the setup, press  one time and select "SAMPLING RATE" with  . Now, choose "MANUAL".
- 5 To go back to the measurements press  three times. Start measuring with  and press  every time you want to record a data.
- 6 Once you are finished measuring stop the Labdisc by pressing  (you will see the instruction "Press SCROLL key to STOP") and press  .

- 1 First, make a sunblock testing-card cutting two windows out of the cardboard and pasting on to it the plastic film. Cover one window with one of the sunblocks and write the SPF in the cardboard surface. Repeat the procedure in next window and label it.



- 2 Place the Labdisc in an open field and record the UV radiation pointing out the sensor directly to the sun.


- 3 Repeat the measurement, covering the sensor with each window of the sunblock testing card. Be sure the sunblock layer on the plastic is facing the sun.
- 4 Measure the UV radiation through the sunglasses, covering the UV Labdisc sensor with two lenses from different sunglasses.
- 5 Once you have finished, stop the Labdisc.

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Experiment

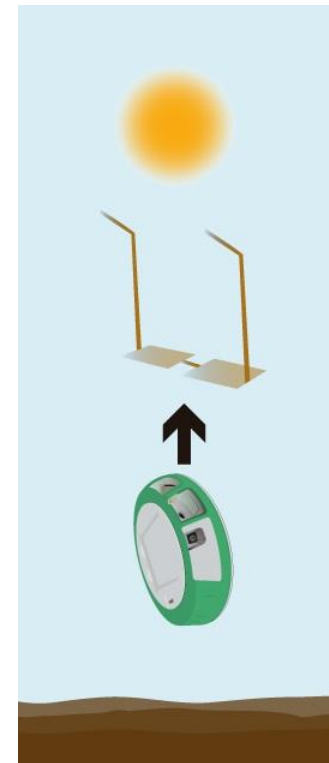





Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Results and analysis




- 1 Connect the Labdisc to the computer using the USB communication cable or via the Bluetooth wireless communication channel.
- 2 On the upper menu press  and select . Choose the last experiment of the list.
- 3 Select a bar graph  from the GlobiLab menu to show the experiment results and label bars using the tool .

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Results and analysis


-  According to the experiment results, do the different solar protection methods work?. Which show better results?
-  Did you find differences between both sunscreens and both sunglasses?
-  How are the results related to your initial hypothesis?

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Results and analysis

The graph below should be similar to the ones the students came up with:

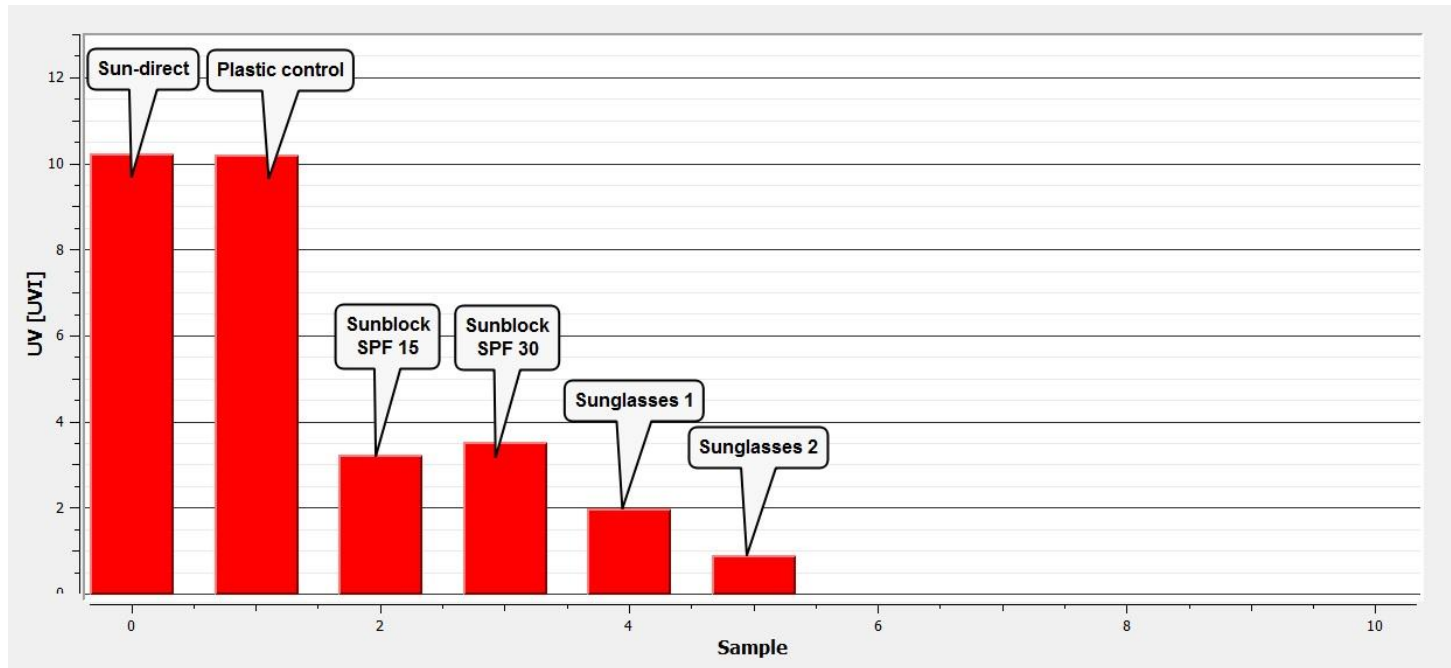

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Conclusion



Why can sunscreen and sunglasses prevent sun damage?

Students should indicate based on the theoretical background that the composition of sunscreen and sunglasses includes chemical substances, which absorb and reflect UV radiation range and other electromagnetic waves, acting as a filters.



Why did the sunblocks and sunglasses show different effectiveness?

Students should infer that differential results were produced by the application of the cream over the plastic. Probably, sunblock is not as effective as on the skin, because the minimum percentage of protection known is around 90%.



Why didn't the sunblocks show a correlation between SPF and UVI measured?

Students should point out the sun protection factor indicates precisely the time of protection offered by the sunscreen. In other words if it takes 20 minutes for your unprotected skin to start turning red, using a SPF 15 prevents burning 15 times longer. However, both types of sunblocks are able to filter out similar percentages of UVB rays (i.e. SPF 15 = 93%; SPF 30 = 97%).

Applied Sciences

UV & Sunblock

Measuring and comparing the level of ultraviolet radiation, through different types of filters.

Activities for further application

? What do you think UV levels depend on? Research and answer the question.

Students should mention some environmental factors related to this phenomenon. It is well known that the ozone layer is the classic guardian against radiation, especially UV. In addition, other factors such as time of the day and time of the year are related to the angle of the sun rays. We can also mention latitude and altitude. For Latitude, at the Equator the UV rays must travel the least distance through the atmosphere. For altitude, high altitude entails less atmosphere to absorb the radiation- and weather conditions.

? UV rays help us to produce vitamin D3, could you propose other benefits offered by UV radiation?

Students could mention the opportunity to use UV radiation as a water purification method. For example, highly efficient mercury lamps transmit a broad range of the EM spectrum, including UV radiation. Pathogen organisms exposed to this then perish.

efectoeducativo

globisens



 **Globisens**
Lab classes with sensors
Labdisc

Digital content
provided by
 efecto educativo
efectoeducativo.com